
DEBLISTERING MACHINE

SIMPATY - F140

Semi-automatic Machine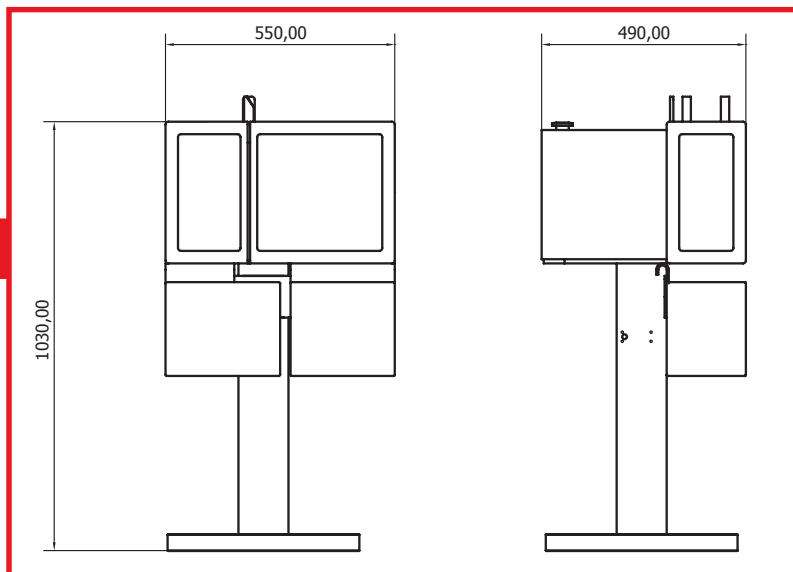

• TECHNICAL DATA •

Max. blister width	mm 120
Max. Blister lenght	mm 200
Standard voltage	230V monophase + E 50 Hz
Dimensions	mm. 620x580x1220h
Weight	Kg. 125
Production	60 blister/minute

• DATI TECNICI •

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- Alu
- PP
- Medical Paper
- Aluminum PE
- PE

APPLICATION:

- Pharmaceutical
- Homeopathic
- Veterinary
- Dental
- Herbal
- Integrators
- Food
- Cosmetic
- and others

THE SEMIAUTOMATIC DEBLISTERING

The semiautomatic deblistering machine SYMPATY F140 distinguishes itself in the pharmaceutical production world for its universality. The thirty-years of O.M.A.R. experience brought about the invention of a unique machine able to deblister all kinds of blister with different formats in plastic material and hard aluminum covering, and ALU-ALU blisters without purchasing any additional tools and with perfect results in terms of product integrity.

UK Sympaty F140 is equipped with all accessories to recover capsules, multi-shaped tablets, soft gel capsules, almonds and so on from blisters with up to 7 pocket rows and a max. width mm 120, thus reducing at zero costs for format parts and their storage. Complete with movable support base in stainless steel and transparent Pet-g bins for the de-blistered product and waste material, Sympaty F140 is supplied with fully interlocked safety guarding with a carter in aluminum and Pet-g that can be open. The recovery bins are fully interlocked to the safety circuit and if removed, or with open carter, the machine automatically stops. The machine can be supplied with electronic piece-counter and blister guide equipment on size for blisters with special configuration such as diagonal placed pockets. The pharmaceutical products are recovered without any damage or contamination thanks to a rotating blade system supplied with Sympaty F140 which cuts the aluminum covering of each single pocket before the product's ejection, in order to help this operation and guarantee the product's integrity.

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